
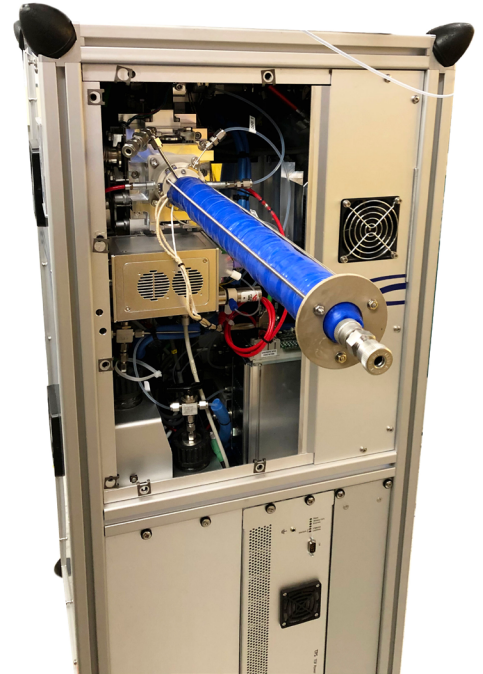


# VIA

## Vaporization Inlet for Aerosols

*Real-time chemical analysis of aerosols with Vocus technology.*


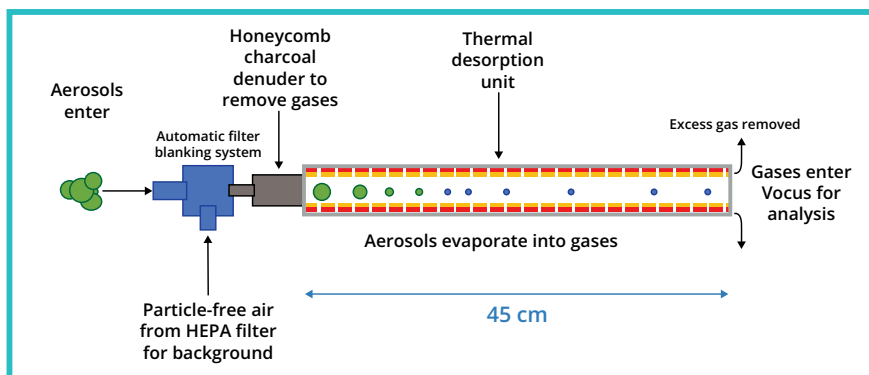


### Applications

- Online identification and quantification of trace-gas- or particle-phase compounds
- Thermal desorption of organic aerosols for analysis by Vocus
- Quantification of organics using  $\text{H}_3\text{O}^+$  PTR-MS
- Detection of HOOM with  $\text{NH}_4^+$
- [egusphere-2023-1146.pdf \(copernicus.org\)](#)
- [acp-23-3707-2023.pdf \(copernicus.org\)](#)

### Advantages


- Lack of aerodynamic lens means small particles  $\text{Dp} < 80 \text{ nm}$  can be measured with the same efficiency as larger particles (no complicated corrections; see below)
- Valve system to enable automated filter blanks and alternating gas- and particle-phase sampling
- Easy to remove and install (< 10 minutes)
- Integrated into Vocus control software

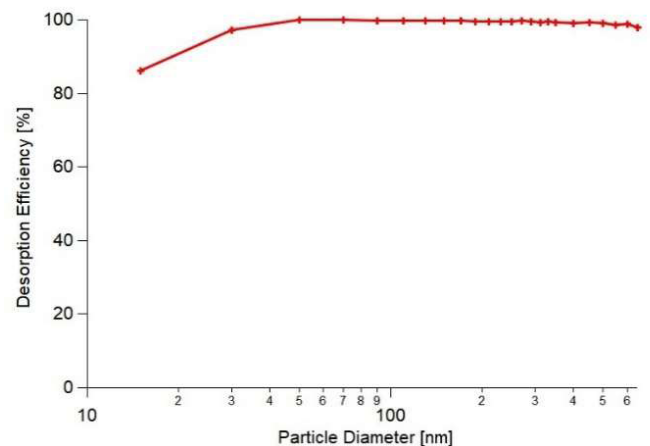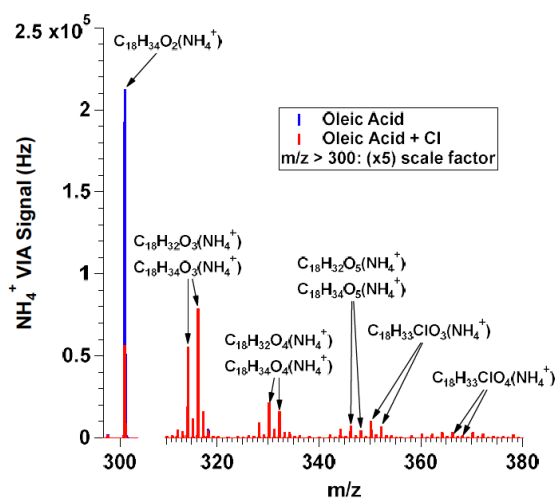

# VIA - Vaporization Inlet for Aerosols

## Performance and Specifications

- ng m<sup>-3</sup> detection limits, without the need for complicated aerodynamic lens due to the high sensitivity of the Vocus PTR-TOF
- Efficient desorption from 25 to 600 nm particle sizes
- Honeycomb charcoal denuder removes > 99% of VOCs while passing through particulates and maintaining sample humidity
- Programmable desorption temperatures from 150 – 250 C, capable of thermogram ramps
- Capable of switchable gas/particle-phase analysis
- Fully integrated into ToFDaq acquisition software or EyeOn control

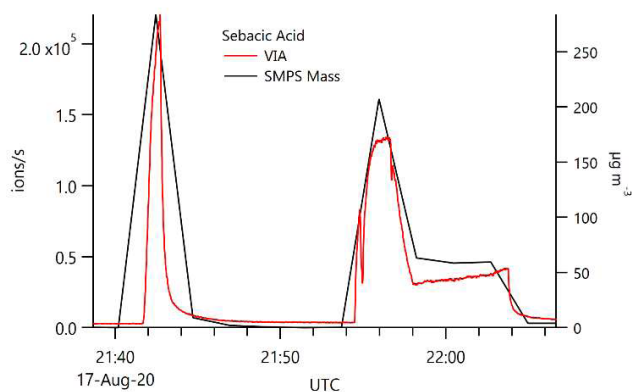

- Left: sebacic acid concentrations next to
- Fast time response to large changes in concentration and has relatively low carryover
- Typical aerosol sensitivities up to ~ 1000 cps/µg using NH<sub>4</sub><sup>+</sup> as the reagent ion

- Automated filter blanking system enables fast and reproducible background determination
- Background periods are recognized and automatically analyzed by Tofware analysis software package
- Below: a time series of azelaic acid with automated filter blanks every 30 minutes

